

Guard columns quick installation instruction

An analytical guard starter kit contains:

- 1 coupler
- 1 column holder (2 parts)
- 5 cartridges

- 1 - Insert the **coupler** at the Kromasil column inlet and finger tighten the coupler onto the column.
- 2 - Insert the **guard cartridge** into the column holder body. Hand tighten the holder lid onto the holder body. Initially hand tighten, then using a pair of 9/16" wrenches tighten a further 1/4 turn to seal the holder lid onto the holder main body (the spanners should grip the holder lid and holder main body).


It is recommended that this seal is not disturbed until the guard cartridge reaches the end of its useful life.

- 3 - Attach the holder to the coupler already mounted on the column inlet and finger tighten.

For latest instruction version and further details, visit:

www.kromasil.com/guards

Holder (cartridge inside) and coupler assembled on column

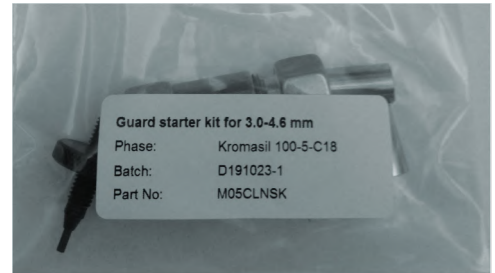

the coupler



the cartridges



the holder

